Characterization of Nanoparticles by Dynamic and Static Light Scattering Technique – [CHP]

(Kutorglo Edith Mawunya, Miroslav Šoóš)

1. Introduction

Light scattering is a powerful non-destructive technique for material characterization, covering size and shape of particles, molecular weight of polymeric chains up to the structure of aggregates formed by aggregation process. The technique is based on the scattering of light by the dispersed particles and depending on the applied setup we refer to the dynamic light scattering (analysis of the scattered light intensity as a function of time) and static light scattering (analysis of the scattering as a function of the scattering angle).



Fig. 1: Cartoon of the Dynamic Light Scattering setup. The key feature is the analysis of time profile of the scattered light intensity measured at a specific angle



Fig. 1: Cartoon of the Static Light Scattering setup. The key feature is the analysis of the scattered light intensity as a function of the detector position. At every angle commonly the average intensity over certain period of time is considered

2. Theory

2.1. Dynamic Light Scattering

In a colloidal dispersion suspended particles undergo Brownian motion. The motion results in fluctuations of the distances between the particles and hence also in fluctuations of the phase relations of the scattered light. Additionally, the number of particles within the scattering volume varies in time. The end result is a fluctuating scattered intensity $I_s(q; t)$. The corresponding measured intensity correlation function is written as:

$$g_{2}(q,\tau) = \frac{\left\langle I_{s}(q,t)I_{s}(q,t+\tau)\right\rangle}{\left\langle \left|I_{s}(q,t)\right|^{2}\right\rangle}$$
(1)

with τ being the time delay. By means of the Siegert relation $g_2(q; \tau)$ can be related to the field correlation function $g_1(q; \tau)$:

$$g_{2}(q,\tau) = 1 + \beta |g_{1}(q,\tau)|^{2}$$
 (2)

with β is so-called intercept. The field correlation function may be used to determine the diffusion coefficient *D* of the scatterers. The field correlation function $g_1(q; \tau)$ can be described by the following relation:

$$\ln[g_1(q,\tau)] = -Dq^2\tau + \frac{c_v}{2}(Dq^2\tau)^2 + \dots$$
(3)

where D is the intensity weighted diffusion coefficient, c_v is the coefficient of variation related to the sample polydispersity. Using the Stokes-Einstein equation and the viscosity of the surrounding medium the radius can be calculated:

$$R_h = \frac{k_B T}{6\pi\mu D} \tag{4}$$

2.2. Static Light Scattering

In static light scattering experiments the intensity of light scattered due to the presence of objects (scatterers) with refractive index different from the surrounding medium is measured as a function of the scattering angle, θ . This angle is directly related to the modulus of the scattering vector $q = 4\pi/\lambda \sin(\theta/2)$, by the wavelength λ . In the simplest case of monodispersed spherical particles the average scattered intensity $I_s(q)$ is proportional to the particle form factor P(q) as follows

$$I(q) = I_0 CNV_p^2 P(q)$$
⁽⁵⁾

where I_0 is the incident light intensity, C is the instrument constant which is a function of the experimental setup and N is the number of particles with a volume V_p . In the case of monodisperse particles with diameter substantially smaller that the laser wavelength P(q) can be calculate analytically as follows:

$$P(q) = \left[\frac{3\left(\sin\left(qR_p\right) - qR_p\cos\left(qR_p\right)\right)}{\left(qR_p\right)^3}\right]^2$$
(6)

with Rp being the particle radius.

Since the size of particles is a priory not known, commonly a Guinier plot is used to extract the particles size. The Guinier Plot allows the determination of the Radius of Gyration (R_g) from the measured scattered intensity $I_s(q)$ as a function of the scattering angle θ . It is used in X-Ray, neutron and static light scattering (SLS). Using a logarithmic plot of the Guinier equation

$$\ln I_{s}(q) = \ln I(0) - R_{g}^{2} \frac{q^{2}}{3}$$
(7)

where R_g can be simple obtained from the slope of the liner fit of the ln $I_s(q)$ vs. $q^2/3$

3. Aims of the work

- Analyze the size of particles synthesized either by emulsion polymerization or precipitation methods.
- Compare the size obtained by dynamic and static light scattering technique
- Investigate the impact of salt concentration on the measured size and stability of obtained particles.

3.1. Pre-lab preparation

Read the laboratory protocol carefully to understand the procedure. Furthermore get familiar with the instrument setup and controlling software.

4. Experimental procedure

4.1. Materials and methods

 Chemicals: Deionized water Sodium Chloride (NaCl), MW: 58.44 g/mol Particle suspension prepared either by emulsion polymerization or precipitation method
 Apparatus: Beakers (50 ml) Pipets (200 μL and 1 mL) Balance

4.2. Work Procedure

- **1.** Prepare a solution of nanoparticle suspension in water. Use 50 mL of deionized water into which you add 200 μ L of original particle suspension.
- **2.** Transfer 2 mL of this diluted suspension of particles to the measurement cuvette and clean its outer part with the paper tissue.
- **3.** Insert cuvette into the device goniometer and select either Dynamic or Static Light Scattering mode. No for should be spplied when inserting the cuvette into the goniometer! It could cause damage of the instrument. Please follow closely the manual or ask the laboratory instructor.
- **4.** Analyze the prepared samples and record the measured values of particle size as well as polydispersity.

5. Evaluation and Report preparation guidelines

This work is supporting the previous step where students prepared particles either by emulsion polymerization or by precipitation. Therefore obtained results will be part of the report for these two works. Protocols including sizes of the prepared particles as well as trends as a function of the condition will be provided not later than 1 week after characterization laboratory described in this document.

6. Precautions and safety information

- (i) **Safety first!** Personal protective equipment (lab coats, gloves, goggles, shoes that cover the entire feet: no slippers or sandals allowed) should be worn at all times when in the lab and when handling all chemicals.
- (ii) information (Source: Sigma Aldrich material safety data sheet) regarding all reagents used in this lab is as follows:

Chemical	Signal word	Symbol
NaCl	No toxicity data available	

Prevention is better than cure!

(iii) Before starting the experiments, ensure all glassware are clean, contamination can alter the results of your hard work.

7. Control questions

- **1.** Which methods for particle characterization do you know?
- **2.** What is the difference between static and dynamic light scattering? How the measurement is done using both techniques?
- **3.** How to calculate hydrodynamic radius and radius of gyration? What is the difference between them?
- 4. What is the impact of the salt on the size of particles?

Good luck!

8. Appendix – Extract from Instruction Manual to the LS Instrument device

8.1. LS Spectrometer software

Version: 6.3 to 6.6

If an older version is installed on your system you might ask LS Instruments for an update.

8.2. Structure of the main program

This section explains in detail the structure and functionality of the LS Spectrometer software provided with the instrument. The graphical user interface is organized into 3 different sections as shown in Fig. 8.1.



Fig. 8.1: Overview of the LS Spectrometer software

The upper section (A) informs about the current status of the measurement and setup: The upper section shows on the left the temperature of the sample. This temperature is based on the temperature reading of a thermometer close to the sample and a calibration which relates the temperature of the sample to the read out of the thermometer. The steady state of this temperature is indicated above with a colored led (grey=unavailable, green=steady state, red=steady state not reached). The two fields below inform about the set values for the sample viscosity and the refractive index of the sample. In the second column, the status of the detector(s) is shown by an indicator (green=on, red=off, grey=unavailable). Below the intensity of the laser is shown, as well as the actual value of the scattering angle and the actual mode of the instrument.

A precise description of the different modes can be found below. In the middle of the upper panel, the actual count rate, its average and standard deviation are shown for both channel separately (if second detector is installed). On the right the timing of the current measurements are indicated. Also the main exit button is located here (Please use only the exit button and not the cross button in the top right corner of the program to exit the)

The middle section (B) is divided into three tabs, *Settings, Measurement* and *DLS Analysis*: This section contains all the software controlled settings of the *LS Spectrometer*. Depending on the hardware configuration and options installed, certain parts are grayed out. The settings are divided into three columns. On the left, measurement settings, in the center the goniometer settings and on the right the laser attenuator settings can be adjusted.

The lower section (C) informs about the initialization and status of the hardware and the software: When the software starts it initializes sequentially all the available options, the corresponding status is indicated by a colored led for each option and details are provided in the status text field on the right. Options which are not available remain grayed-out. A correctly functioning option turns to green after the initialization, while options that fail to initialize are indicated by a red led. When the initialization process is finished, *Ready* is displayed in the text field on the right.

8.3. Measurement settings

8.3.1. Scattering geometry

Here one can select the scattering geometry, which can be either 2D or 3D. In section 3.5 the switching between scattering geometries is explained. Upon the setting change the software will show a small animation illustrating the required corresponding hardware operations. The different options of the selected scattering geometry can be selected in the pull down menu below. For each scattering geometry, there are different correlation types available. The available correlation types depend on the options you have installed and the actual selected scattering geometry.

8.3.2. Scattering geometry and correlation type

1. auto: The auto-correlation is the standard correlation mode. In this mode a single APD is used and the correlation function is calculated auto correlating the corresponding signal.

2. pseudo (optional requiring two APD): In this mode of measurement, the signals of both APDs are correlated. It is only available in the 2D scattering geometry. The pseudo cross mode suppresses after pulsing which has a particularly negative effect on the measurement of small particles. For more information study section 1.

3. 3D Cross (optional requiring 3D option): The 3D cross correlation mode is required for turbid samples to suppress multiple scattering. For more information study section 1 and 2.

4. Mod3D Cross (optional requiring the 3D and the modulation option): Further improvement of the 3D cross-correlation mode. For more information study section 1 and 2.

Auto save data. The "auto save data" option, if checked, allows automated saving of the recorded data while measurements are performed. Click on the folder icon and chose the folder path where you want your data to be saved and a base file name. Consecutive measurements will be saved in

different files having the specified base name followed by running number. The first number of this running number can be specified on the main panel of the measurement settings.

Number of runs per measurement and run duration. The length of a measurement can be defined by two numbers. One is the duration per run and the other is the number of runs per measurement. The total duration of one measurement is thus the number of runs times the duration per run. The splitting of one measurement into several runs has the advantage of allowing to assess the stability of each measurement.

Setup script. The script window is very useful if several measurements in a row should be conducted. A screen shot of the script window is shown in Fig. 4.2. The upper line, "*New Set*", is the line which can be edited. It contains the lowest angle and the highest angle of the series of measurement, the angular step size, the number of measurements and its duration as well as the waiting time in between each run and the temperature. Note that depending on the angular step size the last measurement will be at a lower angle than the highest specified angle. For example: lowest angle 20 degree, highest angle 135 degree and step size 10 degree would lead to an highest angle in this series of 130 degree. Once the parameters are set, you may click on "*Insert line*" to include the line just prepared in the script.

All sets of runs are shown as a list in the middle of the window. If you would like to delete a line click on it and select "*Delete line*". It is not possible to directly edit a line in the main window. On the right panel, the estimated time for the script to be completed is shown.

If this time is unexpectedly large, check the number of runs per measurement as this influences the total script time.

The buttons "Load script" and "Save Script" enable you to save and load scripts. The lower panel in this window shows the file name and location as well as the start number added to the files for auto saved data. Once the script is configured, you may click on the OK button to return to the main window.

Laboratoř přípravy nano a mikromateriálů

I35.0 5.0 1 20 0 24.58 Lower Angle (°) Upper Angle (°) Step # of Measurements Duration (s) Wait (s) Temp (°C) Z 15.0 140.0 5.0 4 40 0 24.58	Delete line	58	24.			weasurements	# of	Step	Angle (°)	Upper	Angle (°)	ower
Lower Angle (°) Upper Angle (°) Step # of Measurements Duration (s) Wait (s) Temp (°C) # 15.0 140.0 5.0 4 40 0 24.58	Delete line			0	20		1	5.0		135.0		30.0
15.0 140.0 5.0 4 40 0 24.58		Toma (%C)	Mait (a)	tion (c)	ontr Du	# of Moscurom	Cton	ade (%)	Unnor A	ala (?)	Louver A	ot #
	Estimated tim	24.58	0	ition (s)	40	4	5.0	igie (*)	140.0	igie (*)	15.0	et#
	01:09:20		-									
	Load Script											
	Save Script											
Name for Auto Saved Data									ved Data	uto Sa	lame for /	File N
									nts\	nstrume	Users\LS I	_ C:\

Fig. 8.2: Overview of scripted window

CONTIN settings. Within a scripted measurement, CONTIN [10] analysis can be performed automatically. If ticked, the settings of the CONTIN analysis can be adapted. The according window is shown in Fig. 8.3. As explained in section 4.1.8 the scattering model, the type of boundary selection, and the number of nodes can be selected. Additionally, the definition of the baseline can be adjusted. This is adjusted in terms of number of channels starting from the highest available channel. For more information on the other parameters please study section 4.1.8.

Scattering Mod	del	C et Dev	
R Min	R Max	Set Bo	unus manualiy
1E+0 nm	1E+3 nm		
Nodes 150 🗘			
Number of Ch in the Baseline	annels		
10 🗘	Save	Contin In/Output	t FIles

Fig. 8.3: Overview of scripted CONTIN analysis settings

Channel scaling method. When performing measurements on the same sample but at different angles, the measured correlation function shifts along the lag time abscissa. This can cause problems for the cumulant analysis fit. To avoid this issue you can apply lag time scaling. Upon an angle change, the software will try to guess an optimal channel interval for the new angle depending on the chosen channel scaling method.

- 1. **No scaling**: No scaling is applied. The maximum lag time for the fit is constant for all measurements.
- 2. **q squared**: The adjustment of the maximum lag time follows from that of the previous angle by the theoretical q2 law. Details on setting the maximum lag time are described in section 4.1.7.
- 3. **Decay factor**: The adjustment of the maximum lag time is based on the decay of the g2 function. The maximum lag time is thus set at the lag time where the correlation function decreases from the minimum set lag time by a given percentage which can be fixed by the user. Details on setting the decay factor are described in section 4.1.7.

Total counts threshold. This value is used for the cumulant analysis performed during the measurement. A minimum number of counts is required to perform a cumulant analysis with a meaningful result. The total counts threshold fixes this threshold.

Manual input temperature. If no software controlled thermostat is attached to the system, this input option appears in the program. Use it to manually set the temperature of the sample.

8.3.3. Goniometer settings

The goniometer settings control the movement of the main goniometer and sample goniometer (optional). For the main goniometer chose a scattering angle and move the arm by clicking "Go". The sample goniometer has a drop-down menu with four entries.

- 4. **Off**: Switches the Sample Goniometer off.
- 5. **Square Cell**: For usage of square cells. The angle of the square cell is adjusted according to the angle of the main goniometer following the θ 2 θ scheme [6].
- 6. **Non-Ergodic DLS**: Sample rotation speed is optimized for best results of a non-ergodic sample in a DLS measurement.
- 7. **Non-Ergodic SLS**: Sample rotation speed is optimized for best results of a non-ergodic sample in a SLS measurement.

8.3.4. Sample properties

This section describes the relevant physical properties of the solvent used for the sample. The solvent can be chosen from a list of standard solvents. This list contains: Water, Methanol, Ethanol, Toluene, Glycerol.

For any other solvent the viscosity and the refractive index can be defined by the user. For this purpose either choose "other" or "load file" from the list. The option "other" allows the user to directly enter the viscosity and the refractive index. The option "load file", enables specification of an arbitrary temperature dependence for the solvent refractive index and viscosity by means

of a text file. The corresponding file format is illustrated in the following example. The sign tab corresponds to a *tab* and *eol* corresponds to a line break:

Temperature tab Refractive Index tab Viscosity (cP) eol eol

280 tab tab 1.32 tab tab 0.8 eol 285 tab tab 1.33 tab tab 0.89 eol 290 tab tab 1.34 tab tab 0.78 eol 295 tab tab 1.53 tab tab 0.76 eol 300 tab tab 1.23 tab tab 0.75 eol

This text can be saved in a normal text file. Respecting the precise syntax is important for a correct read out. Several advanced text editors, such as for example Notepad++, allow the user to display tabs and line breaks, this greatly facilitates the creation of the file above.

8.3.5. Thermostat settings

The temperature of the sample can be controlled via an external thermostat. The temperature of the sample can be set here as well as the accuracy of a temperature. The options "*low accuracy*" and "*high accuracy*" define when the temperature will be considered sufficiently stable for a measurement. Higher accuracy of course imply larger equilibration times.

8.3.6. Laser attenuator (+Detector power)

This panel controls the automated laser intensity attenuator (optional) as well as the detector power state. The automated intensity attenuator can be controlled in three different ways:

- Auto: In the auto mode the incident laser intensity is set automatically to yield an optimal scattered count rate based on the best range of the detector. Detectors are automatically powered on and off. Before and after a measurement, the beam path into the scattering cell is automatically closed.
- 2. Set Scattered Intensity: This option allows the user to perform an experiment yielding a defined scattered count rate in the experiment. In the field "Target Scattered Intensity (kHz)" you can define the desired scattered count rate. Upon measurement start, the software will find the right laser attenuation. You can also define the precision by which the software should try to obtain the desired scattered count rate. The requested precision should be chosen depending on the sample. As a general rule monodisperse samples will allow the use of higher precision settings, whereas for more polydisperse systems higher precision settings might cause the automatic attenuation algorithm to fail. In such cases the user should choose a lower precision setting. Detectors are powered on whenever they are needed and switched off in the standby. The laser is fully attenuated whenever it not used.
- 3. Set Laser Intensity: This option gives the user full manual control over the laser attenuation. You can either type in a value in terms of percentage of the full laser power or use the slider on a logarithmic scale and then click"Go". If this option is used, detectors have to be powered on manually by using the tick box right next to "Detector power". Make sure that you tick this button only if you can't damage the detectors.

If the automated intensity attenuator is not installed you switch the detectors on and off via the tick box.

8.3.7. Focal length compensation

For the auto focus system (optional) you need to enter the type of scattering cell used in the measurement. For square cells only the optical path length needs to be entered. After a change, the button "*adjust*" needs to be clicked. Starting a measurement will adjust the focal system according to the current settings.

8.3.8. Measurement

The measurement panel (Fig. 8.4) shows all relevant information during a measurement. After defining the parameters in the settings panel, click on the button *"Start Measurement"*. In case you have chosen a scripted measurement this button is replaced by the button *"Start Script"*. The buttons to the right allow you to stop or pause a measurement. The buttons below enable you to save recorded data, if not already saved automatically and to load previous measurements. A summary of the measurement conditions which will be also saved in the file is presented below the control buttons.



Fig. 8.4: Overview of the software during a measurement

Count rate trace: On the upper left side are the count rate traces for both (or one) detector. On the right side of this graph the distribution of the count rate is shown. This latter graph is useful for assessing the quality of a measurement. If, for example, a dust particle enters the scattering volume the distribution will clearly indicate this by displaying a long tail towards large intensities.

Correlation function: The graph displays the measured correlation function. From the drop down menu you may chose between logarithmic and linear scaling of this plot. Scaling of the axis can either be performed by the autoscale button or by clicking on the numeric values of the axis. When, during a measurement, the total counts threshold, as specified on the settings panel, is reached additional black, blue and red line appear on the graph. Those lines define the lowest and highest channels defining the correlation function subset taken into account in the cumulant analysis fit. The red line is used for the lower limit and the black and blue one is used for the upper limit. Depending on the channel scaling method you may either move the vertical or the horizontal line, which is indicated by its dark blue color. The limits of the cumulant analysis can be alternatively defined by the numeric fields below the figure. This can be performed in terms of channels of the correlator or in time steps.

If the scripted CONTIN analysis is performed, an additional light blue line appears in the plot indicating the limit of the baseline the analysis.

Measurement trend / angle plots /time trends (from vs 6.6): The lower figure on the right side can show the angular trends of the intercept as used for SLS. The time trend of the hydrodynamic radius and the intensity can for example be used to observe aggregation and the panel of the 2nd order radius and the intercept as a function of the measuring time can be used for determining the quality of the measurement. This panel also indicates the number of total counts and information of the cumulant analysis as the first, second and third order of the radius and its width.

The appearance of the graphs can be edited by a right click on the figure.

8.3.9. DLS Analysis

The DLS Analysis offers you cumulant and CONTIN [10] analysis of single DLS measurements (Fig. 8.6). A theoretical overview of those analysis are given in the Appendix A.

The panel is divided into two sections. On the left are the entries of the analysis as the plot of the data, the limits and settings of the cumulant and the CONTIN analysis. On the right part, the results are presented. As in the plot of the correlation function on the measurement panel, the plot of the measured g² function is shown. The limits for the cumulant analysis can be modified by the blue and the red bar in the figure by simple drag or via numeric values in the panel below. Additionally, the resulting simulations (CONTIN and cumulant) can be shown in the figure. In the case of the CONTIN analysis the simulation appears only after it was calculated by clicking on the button CONTIN analysis (on the bottom right side).



Fig. 8.5: View of the time trends

Parameters for the CONTIN analysis can be set in the CONTIN Settings box. As scattering model either a RGD sphere or point scatter model can be chosen. The latter assumes the scatterers to be significantly smaller in size than the incident laser wavelength while the former relaxes this assumption. The number of nodes (discrete radii) calculated can be specified, too. This number defines the grid of the calculation. The more points are taken for the grid the higher is the resolution of the result. Finer grids result, however, in longer calculation times. It is also important to choose the grid size accordingly to the quality of the measurement. If not ticked, the boundaries of the radius are defined automatically according to cumulant analysis radius. By ticking "Set Bounds Manually" you might define the smallest and the largest radius of the simulation manually.

On the right top part, the results of the cumulant analysis are shown which can be saved with the "Save Fit" button. On the lower part of this panel, the CONTIN simulation results are shown either as a graphical distribution or as a list of peaks within the distribution.



Fig. 8.6: Overview of the Software for DLS analysis

8.4. Structure of the saved files

Several files can be produced as result of either a measurement or an analysis.

8.4.1. Data files

The output format of all files is pure text, which can be imported into your preferred datatreatment program. A single measurement (here for example made in Pseudo cross-correlation) produces an output as shown below, whereas *tab* corresponds to a tab and *eol* corresponds to a line break:

10/29/2013 tab 12:01 PM eo/ Pseudo Cross Correlation eo/ Scattering angle: tab 15.0 eo/ Duration (s):tab 40 eo/ Wavelength (nm): tab 632.8 eo/Refractive index: tab 1.330 eo/ Viscosity (mPas):tab 0.896 eo/ Temperature (K): tab 297.9 eo/ Laser intensity (mW): tab 0.029722 eo/ Average Count rate A (kHz): tab 334.5 eo/ Average Count rate B (kHz): tab 294.0 eo/ Intercept: tab 0.9316 eo/

Laboratoř přípravy nano a mikromateriálů

```
Cumulant 1st tab 295.04 eol
Cumulant2ndtab303.97tab80.62eol
Cumulant 3rd tab 307.54 tab 117.01 eol eol
Lagtime (s)g2-1eol
1.250000E-8tab0.917454eol
2.500000E-8tab0.928283eol
3.750000E-8tab0.935385eol
...
```

```
••
```

```
Count Rate History (KHz) CR CHA / CR CHB eol
0.000000tab 194.520950tab 184.640884eol
0.104858tab 459.136963tab 411.806107eol
0.209715tab 651.416779tab 559.101105 eol
0.314573tab 204.992294tab 183.868408 eol
```

The first 14 lines are used to summarize the experimental conditions as well as the result of the cumulant analysis. Below the summary values of the $g^2(t) - 1$ function are printed in a tab separated format. At the end of the $g^2(t) - 1$ function an under sampled history of the count rate is shown. The first few lines of such history are also printed above for the sake of illustration. If a script was performed, an additional file is saved. While each separate measurement in a script is saved with a running number, the summary file does not have a running number. Its structure is shown below:

```
Scattering angle Mean CR * sin(angle) / Laser Intensity (kHz/mW) g2(t=0)-1 CR CHA (kHz) CR
CHB (kHz) Temperature (K) Laser intensity (mW) eol
15.004272 tab 2735.121749 tab 0.931612 tab 334.522076 tab 293.984703 tab 297.860000 tab
0.029684 eol
15.004272 tab 3363.434631 tab 1.007546 tab 413.088534 tab 360.682696 tab 297.890000 tab
0.029711 eol
19.997803 tab 2650.292570 tab 0.863937 tab 409.287551 tab 358.341338 tab 297.890000 tab
0.049417 eol
```

```
····
```

It contains thus the scattering angle, the count rate corrected by the scattering angle and the intensity, the $g^2 - 1$ function, count rates for both channels, the temperature and the laser intensity. Note: Scaling of the intensity by measured laser intensity is only performed in case of an installed automatic intensity attenuator (optional).

8.4.2. Analysis files

The file created by the button "Save fit" in the cumulant analysis section will output the same file structure as shown above for a single measurement.

The button "Save Fit" in the CONTIN section outputs a ".fit" file which is also a pure text based

Laboratoř přípravy nano a mikromateriálů

file containing the calculated size distribution. Its structure is described below by an example recorded in Pseudo cross-correlation. (*tab* corresponds to a *tab* and *eol* corresponds to a line break.)

10/29/2013 tab 12:03 PM eol Pseudo Cross Correlation eol Scattering angle: tab 20.0 eol Duration(s): tab 40 eol Wavelength (nm): tab 632.8 eol Refractive index: tab 1.330 eol Viscosity(mPas):tab 0.896eol Temperature (K): tab 297.9 eol Laser intensity (mW): tab 0.049960 eol Average Count rate A (kHz): tab 409.3 eol Average Count rate B (kHz): tab 358.3 eol Intercept: tab 0.8639 eol Cumulant 1st tab 338.67 eol Cumulant2ndtab316.84tab120.52eol Cumulant 3rd tab 317.35 tab 116.95 eol eol radius(nm)rel.intensityeol 31.700000 tab 0.000000 eol 32.700000 tab 0.000000 eol 33.700000 tab 0.000000 eol

•••